



## Broken Glass Disposal

### 1. Purpose:

---

To provide step by step guidance on how to handle broken glass.

### 2. Prerequisites:

---

You are an authorized user of DB440 and are either included in your PI's permit, or you possess a CL2 permit for DB440.

### 3. Responsibilities:

---

It is the responsibility of all faculty, staff and students to follow the procedures described in this SOP.

### 4. Personal Protection Equipment (PPE):

---



### 5. Procedure:

---

#### Broken Glass

- 1) Wear PPE as described above. Use a brush and dust pan (in Spill kit) or tongs, etc. to pick up broken glass pieces.
  - i) **DO NOT** pick up broken glass with your hands.

- 2) Dispose chemically contaminated broken glass in the chemical waste pail. Dispose biologically contaminated broken glass in the Bio waste pail as per the [“Biohazard Waste Disposal” SOP](#).
- 3) Remove and discard your gloves. Wash your hands with soap and water.